

High Throughput Solid Phase Extraction, Cleanup and Concentration of Organochlorine Pesticides in Waste Water Using EPA Method 608.3

Ruud Addink and Tom Hall
Fluid Management Systems
Billerica, MA



- Organochlorine pesticides are man-made organic chemicals with a history of wide spread use around the world
- Persistent to degradation, found in drinking water supplies and sediments
- Some have been added to Stockholm Convention (2009-2015)
- EPA method 608.3 - Waste Water



Health risks OCPs

- Organochlorines have a wide range of both acute and chronic health effects
- Cancer
- Neurological effects
- Birth defects
- Many OCPs are also suspected endocrine disruptors



Analysis of OCPs

- Many labs analyze drinking and waste water samples
- Liquid-Liquid Extraction (LLE) or Solid Phase Extraction (SPE) can be used
- In both cases organics are transferred from water sample to an organic solvent
- With SPE compounds are first deposited on cartridge or disk, then eluted



Comparison of LLE/CLE vs. Semi-Automated SPE Methods (1)

LLE/CLE

Open to laboratory background

Uses >360mls solvent

Shaking / Continuous process

Forms emulsions requiring centrifuging

Little Selectivity

Requires water removal

Semi-Automated SPE

Closed system

Uses <60mls solvent

Filtration process

No emulsions formed

Wide Selectivity (adsorbent)

In-line water removal



Comparison of LLE/CLE vs. Semi-Automated SPE Methods (2)

LLE/CLE

No Separation of waste

More volume to evaporate

Massive solvent emission

CLE uses a lot of power

Requires lots of solvent for cleaning

Semi-Automated SPE

Separates Aqueous and Organic Waste

<60mls solvent to evaporate

6 times less solvent emission

Easily Capture Solvent

Lower solvent costs

Lower Disposal Costs

Reduced Solvent Usage



Objective for Semi Automation

- Use as many features from Automated systems and implement them into a Semi automated platform
- Develop as many SPE procedures for the testing lab using a single extraction platform.
- Minimize manual steps to lessen error and maximize limited man hours



Goal (1)

- **Easy to Operate**

- No Computers or Electronics to fail or maintain

- **Semi - Automated**

- Hyphenates the entire Solid Phase Extraction Process – Extraction, Bottle Rinse, Inline Drying and Optional Direct to GC Vial Concentration

- **Fast**

- The fastest sample processing available for SPE
- Run up to twelve samples simultaneously
- Vacuum for fast loading of large volume samples
- Unattended Sample loading walkaway time

- **Closed system**

- Eliminate potential outside contamination




Goal (2)

- **Efficient**


- Uses all SPE cartridge sizes
- Dedicated manifold for cartridge conditioning and sample loading
- Dedicated manifold for extraction and extracts
- Separates Organic from Aqueous waste
- Vacuum cartridge drying, Nitrogen cartridge drying or combined
- Automated Bottle Rinse and Elution
- Inline Extract Drying
- Small number of components to clean




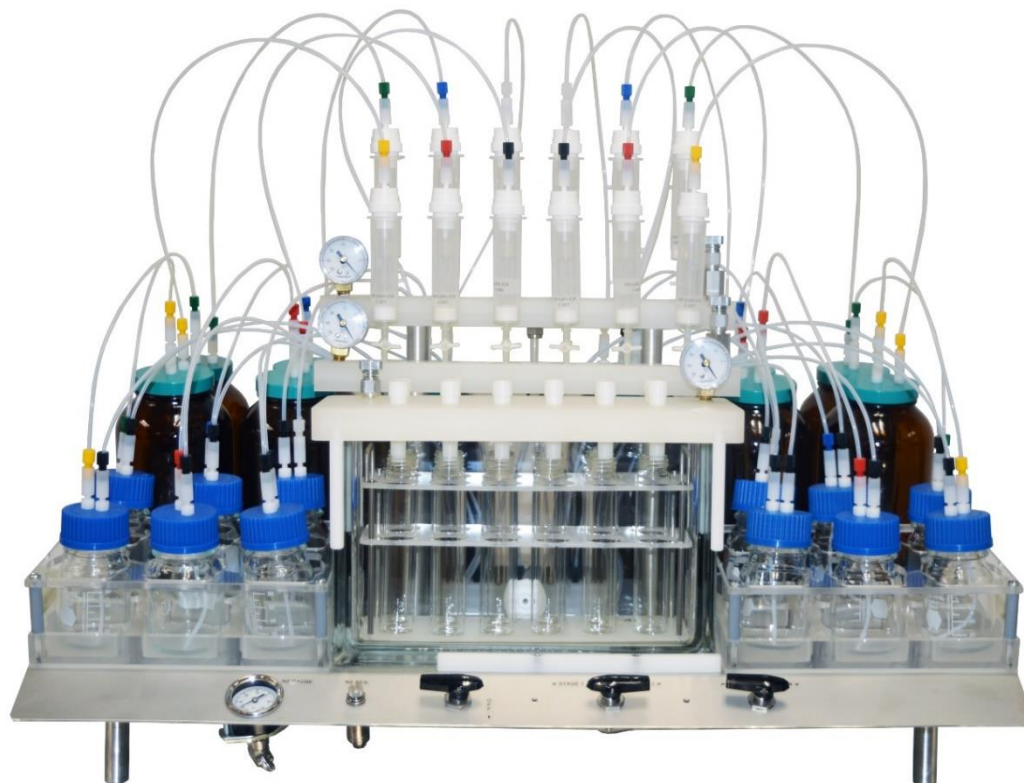


EZSpe 12 samples

System Layout


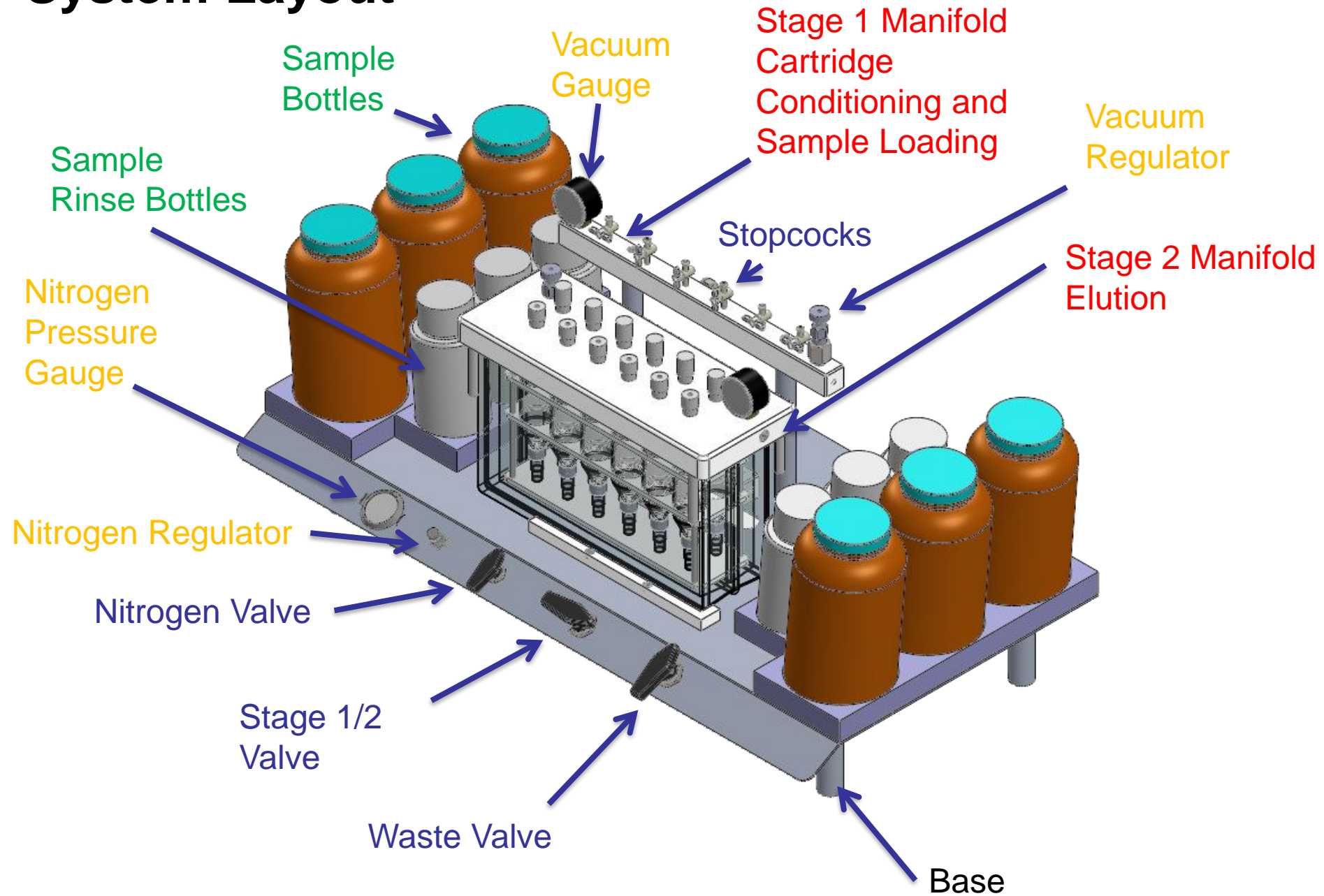


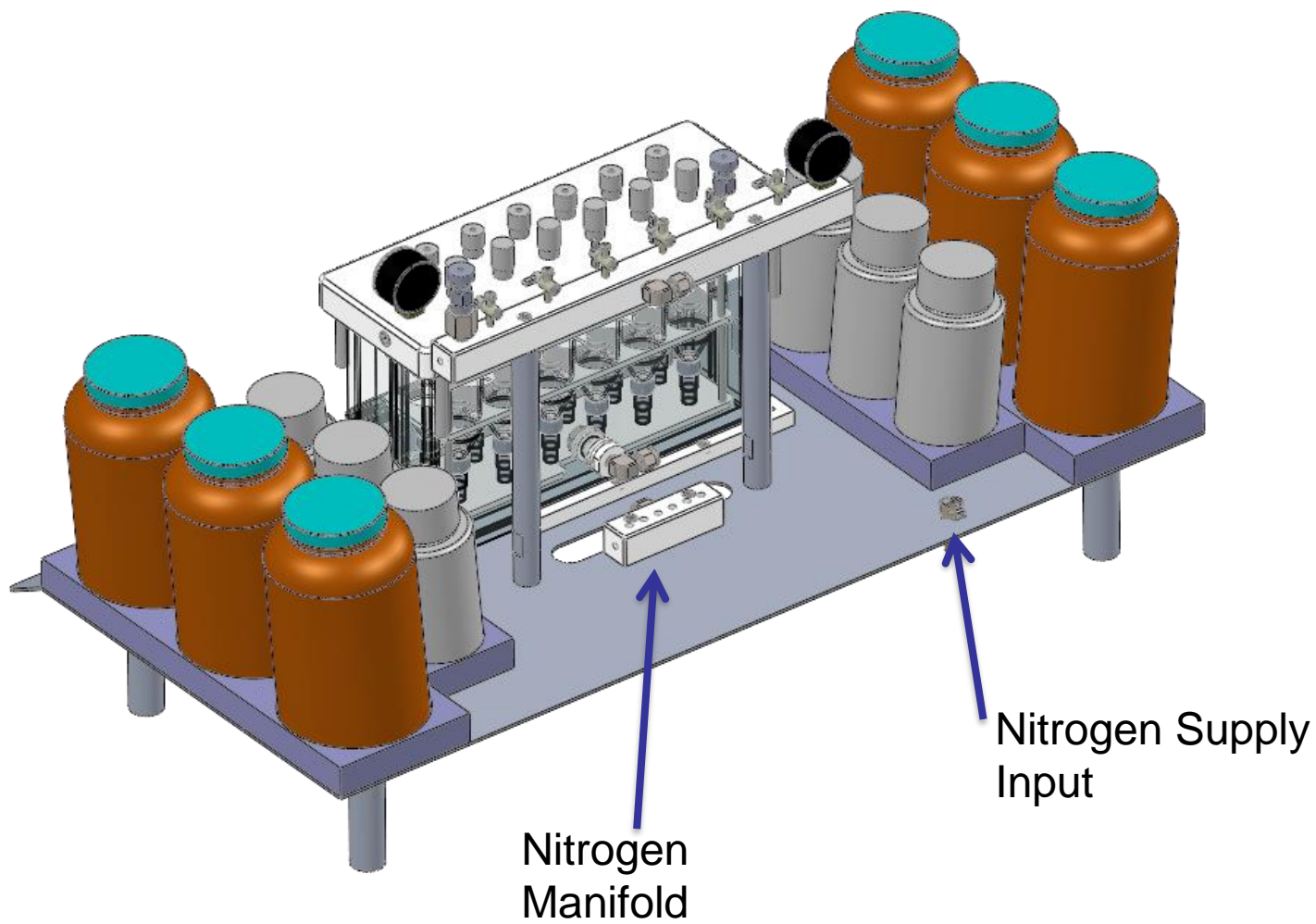


Control Valve Layout

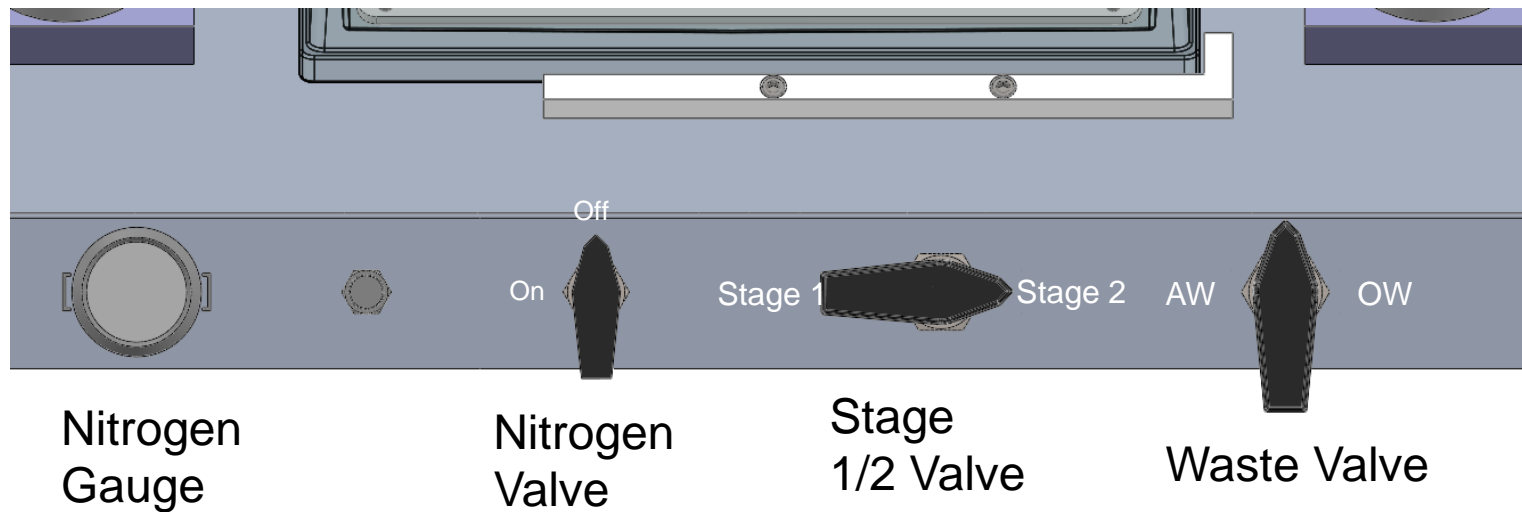


Cartridge Conditioning (Stage 1, Organic Waste)

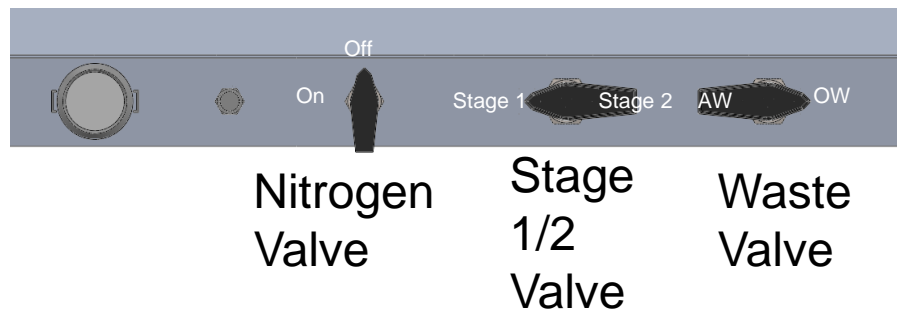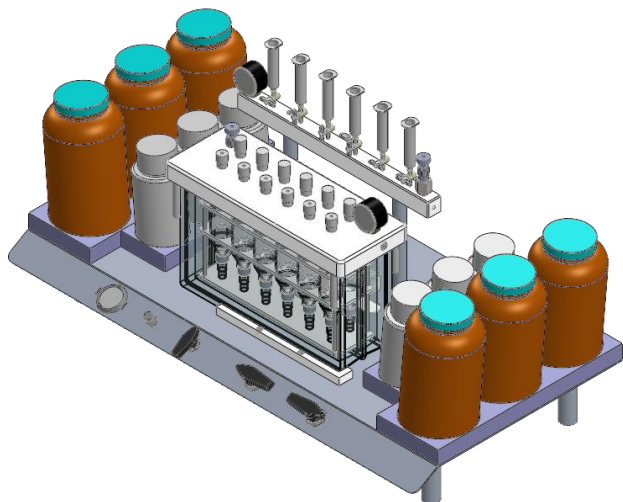


Flow
Path

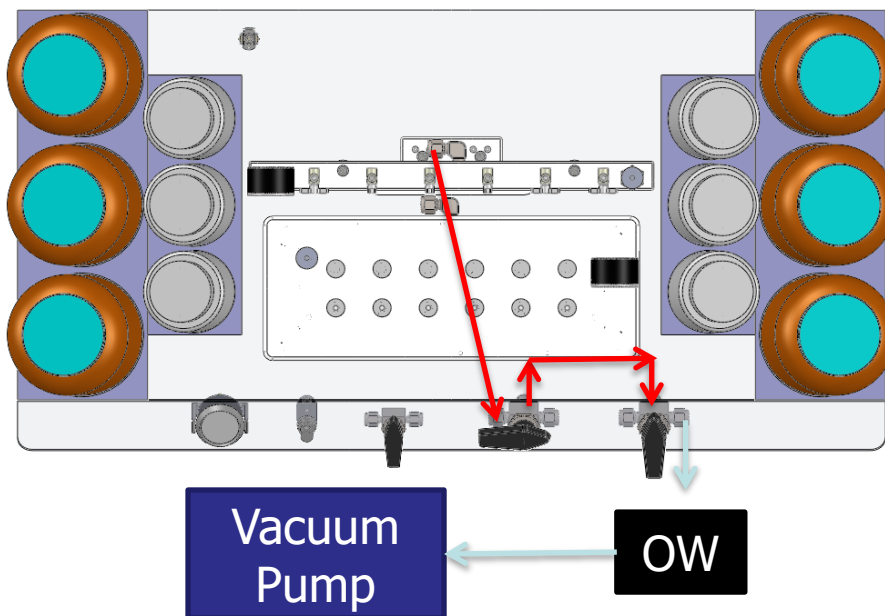

Sample Loading (Stage 1, Aqueous Waste)

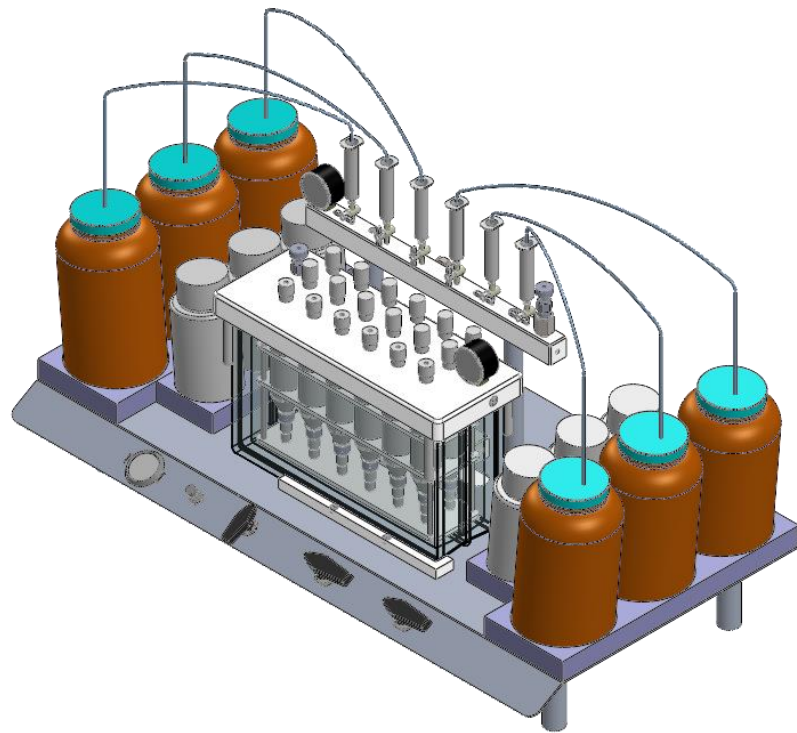


Flow
Path

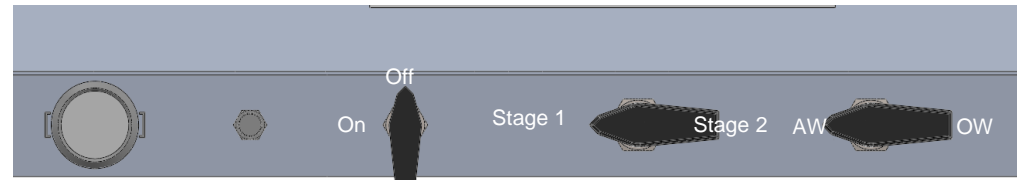

Nitrogen
Valve

Stage
1/2
Valve


Waste
Valve

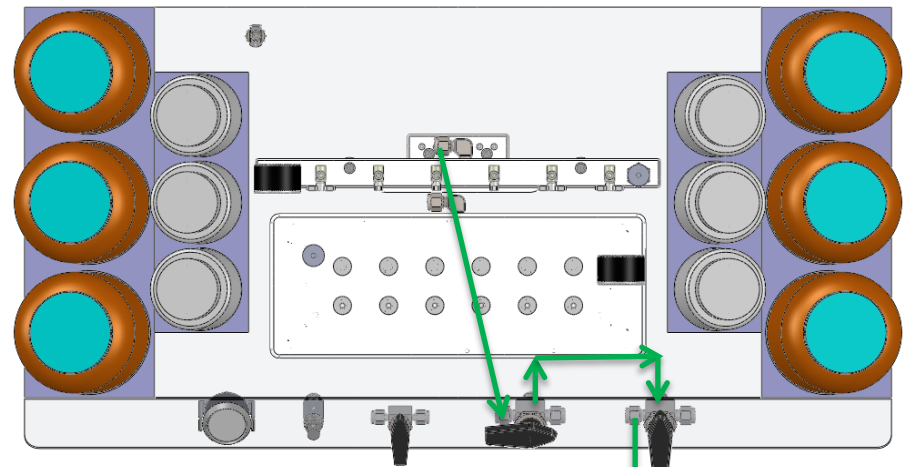


Vacuum
Pump

AW

Cartridge Drying- Nitrogen/Vacuum

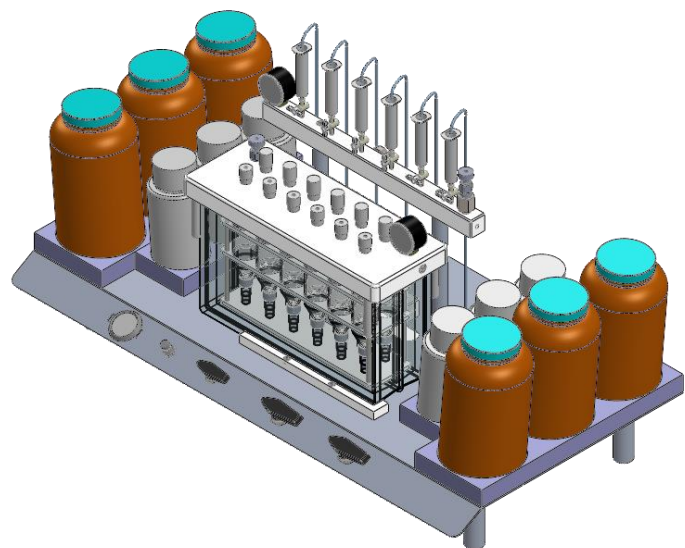


Flow
Path

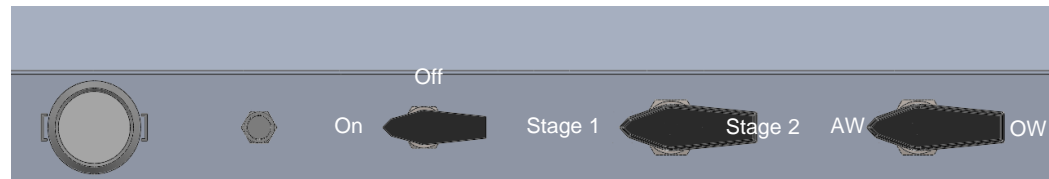


Nitrogen
Valve

Stage
1/2
Valve


Waste
Valve

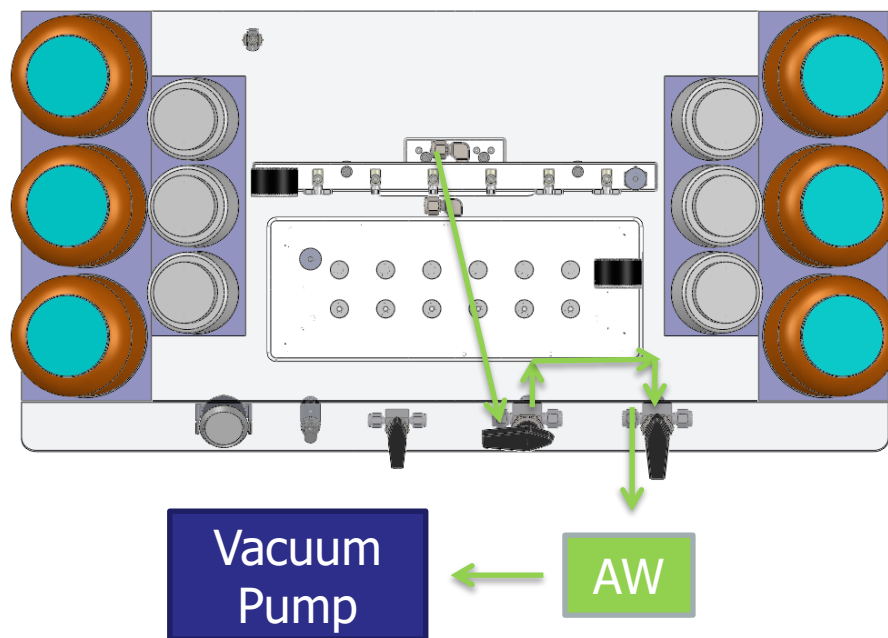

Sample Elution (Stage 2)

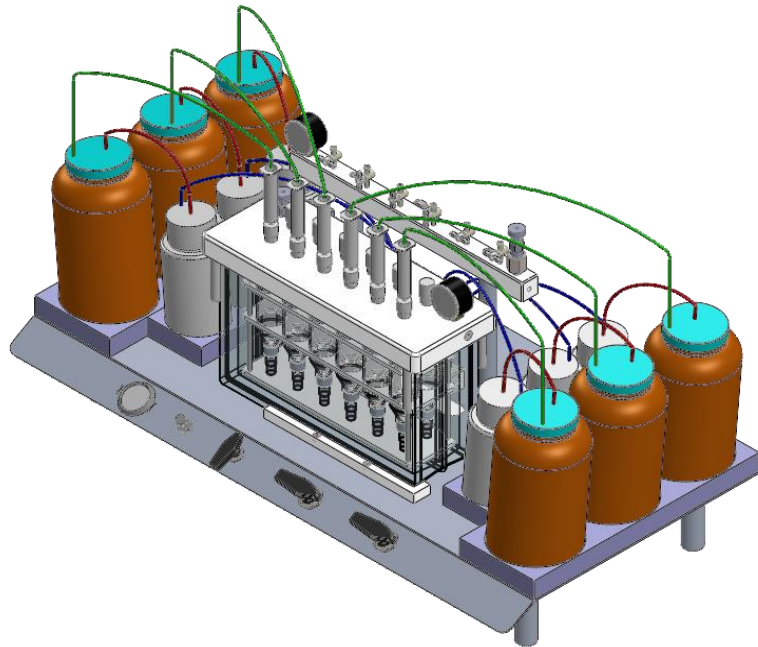


Flow
Path

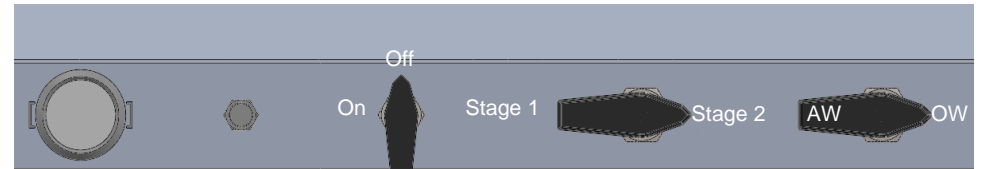

Nitrogen
Valve

Stage
1/2
Valve

Waste
Valve

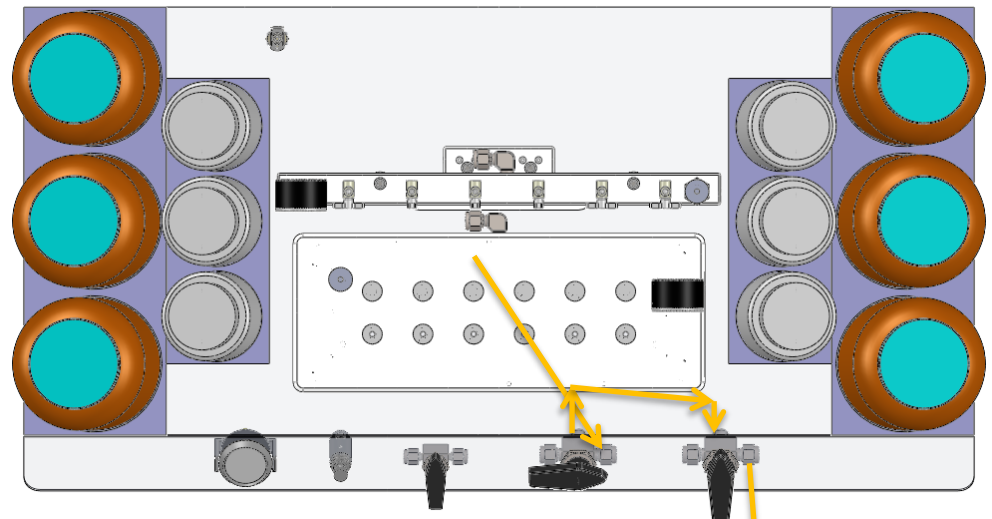

Vacuum
Pump

OW

Procedure (1)

- 12 samples (1L water each) are prepared and acidified with 6 N HCl till $\text{pH} < 2$
- Spike with relevant standards
- Put sample bottles in place and fill automated rinse bottles with 40 mL 10% acetone/hexane
- C-18 cartridges are installed in each of the 12 positions.



Procedure (2)

Stage 1:

- Vacuum is turned on
- C-18 cartridges conditioned with 10 mL acetone (2 min soak), and 20 mL water
- Samples are loaded across cartridges under vacuum (~ 12 inch Hg)
- C-18 cartridges are dried under vacuum for 10 min and then removed
- Hydromatrix cartridges on top of Florisil cartridges conditioned with 20 mL 10% acetone/hexane



Procedure (3)

Stage 2:

- C-18 cartridges on top of Hydromatrix/Florisil cartridges assembly on top of Stage 2 manifold
- Sample bottles are automatically rinsed from the rinse bottles with 40 mL 10% acetone/hexane
- Sample bottles rinses loaded across all three cartridges and collected for analysis (Direct-to-Vial)



Direct to Vial Concentration



SuperVap Concentrator

- 12 position
- 50mL vessel



Concentration Features

- 6 (250ml) and 12 (50ml) position models for extractions.
- Dry bath heating element
- Independent secondary heater for extract nipple (Can be disabled).
- Sensor controlled
- Savable temperature log capability.




Evaporation

- Pre-heat temp: 55 °C
- Pre-heat time: 15 min
- Heat in Sensor mode at 55 °C under nitrogen (7-10 psi)
- Direct to GC Vial Vessel Reduce to 1 mL
- Samples are now ready for analysis (ECD)



608.3 Results

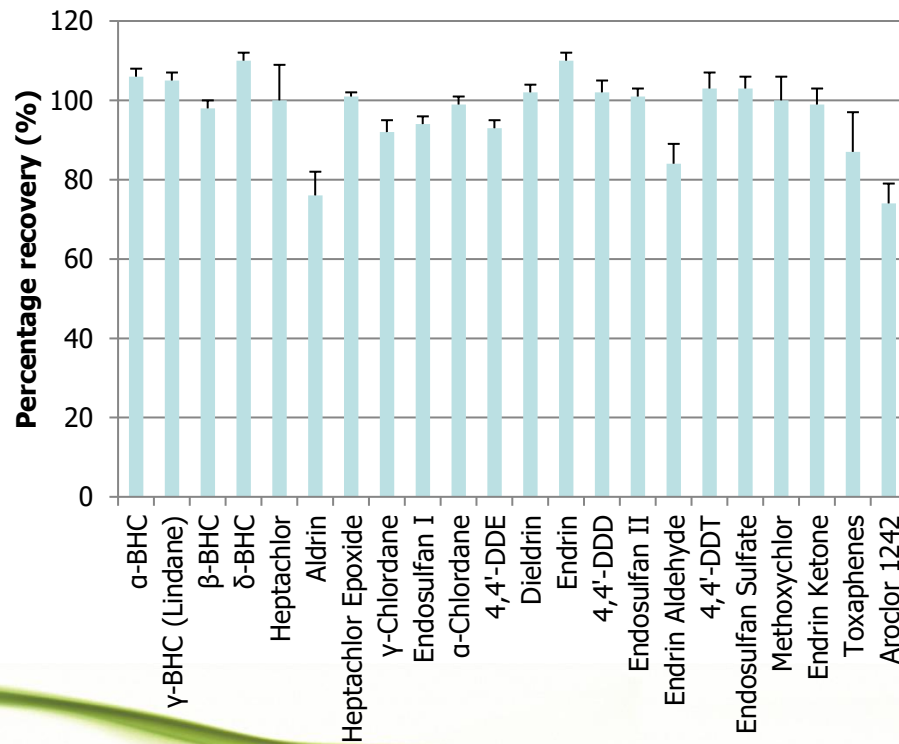

Semi-Automated SPE in Summary

- EZSpe and SuperVap systems are easy to use and install
 - Complete Water Sample Prep Workflow
- Low cost, High throughput, Low maintenance solution
- EZSpe Extractions and Concentration is a very green technique
 - Reduces Solvent Use
 - Reduces Solvent Disposal Costs
 - Reduces Solvent emissions



Semi-Automated Summary

- Semi-automated SPE and SuperVap systems deliver consistent, reproducible results
- Handles a wide range of Sample sizes and matrix types
- Uses all SPE Cartridge and Column sizes
- Comply with existing methods that require vacuum, positive pressure and precise delivery of sample and solvents



Questions

Contact us at:

raddink@fms-inc.com
(Ruud Addink)

thall@fms-inc.com
(Tom Hall)

